**NATURE &
DECOUVERTES**

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BINOCULAR MICROSCOPE

LUPA BINOCULAR

BINOCULAIRE MICROSCOOP

LUPA BINOCULAR

Réf. 42006910

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PLEASE CAREFULLY READ THIS MANUAL AND KEEP IT IN A SAFE PLACE

WARNING

- Non-rechargeable batteries are not to be recharged.
- For electric toys using rechargeable batteries, the batteries should be charged under adult supervision.
- Different types of batteries or new and used batteries are not to be mixed.
- Batteries are to be inserted with the correct polarity (+ and -).
- Exhausted batteries are to be removed from the toy.
- The supply terminals are not to be short-circuited.

DESCRIPTION

This Mini stereo plastic microscope has sharp stereo erect image over a wide field view. Widely used in electric factories, laboratories, schools and families.

TECHNOLOGY DATA

- Upright binocular view head, the interpupillary distance is between 55-75mm.
- The working distance of 2X magnification of objective 70mm.
- Built-in incident light 3V/0.25W.
- Power 2 units of AA batteries (not included).
- Plastic body and easy to carry.

HOW TO USE

- The room temperature for the microscope is between -5°C + 40°C .
- Inserting and removing batteries: remove the battery case cover using a Phillips screwdriver, turning anticlockwise. Then place 2 AA batteries in the case, carefully observing the +/- polarity of each battery. Finally, once all the batteries have been inserted, re-fasten the battery compartment cover using a Phillips screwdriver, turning clockwise.
- Touch switch position "ON", the microscope light will turn on. Place the specimen in the center of the working plate and fix it with stage clips. Loosen fastening knob and move the head up and down until the specimen within the working distance then fasten it. Turning focusing knob slowly until there is sharp image through right eyepiece. Then observe left eyepiece, if the image not clear enough, then adjust diopter tube, Rotating the prism house to suite your interpupillary distance. After the above, a crisp erect image appeared.
- If needed, do not forget to turn-off the light when replacing LED lamp. Take off the lamp house screw and replace the lamp with new LED lamp.
- **Note:** Objective must be on the working stage when using.

MAINTANCE AND REPAIRE

- Put the microscope in a cool and dry place.
- Keep the microscope clean, not forget to put some grease to driving system from time to time.
- All the optical components are adjusted and could not be taken apart by yourself. When the lens and other optical parts covered with dust, you should wipe with volatile solvents or abrasive cleaners.
- Deliver to a professional shop for repair when not in use.

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